


# **Biomaterials Engineering**

M3D1 11.15.11

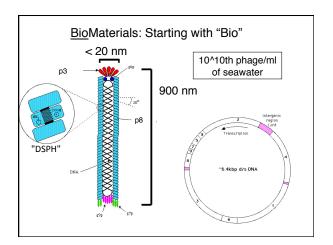

# Phage life cycle 1. Infection 2. Amplification 3. Morphogenesis + release of progeny phage • new M13 phage particles within 10' post-infection • 1000 phage/cell within the first hour

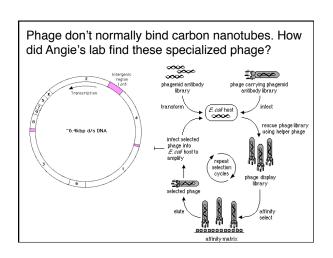

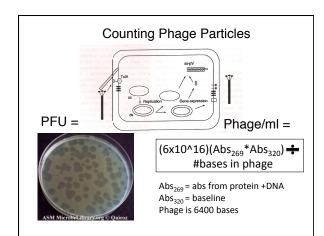


# Parts 1 + 2 of your experiment today


### Phage purification

- · Harvest phage infected cells. Phage in SUPERNATANT!
- Concentrate by precipitation (+PEG/salt)
- · Pellet may not be visible...have faith.
- · Pool phage with "super-group"

## Measurement of phage concentration

- · Use quartz cuvettes
- Dilute 1:10 in TBS and measure Absorbance
- Calculate volume needed for 4x10<sup>13</sup> phage (want this volume btw 0.5 ml and 5 ml)

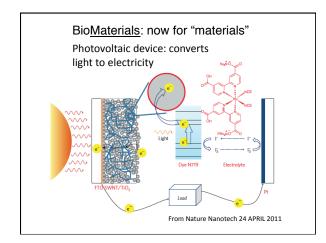

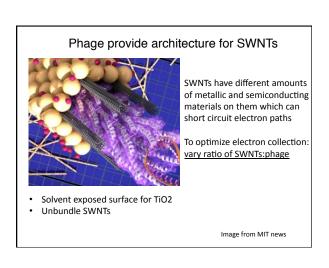

# Part 3: reacting phage with SWNTs

Ratio SWNTs:phage	Super-Team
1:1 (40 ug:4*10^13)	Red, Green, Purple
2.5:1 (100 ug:4*10^13)	Orange, Blue
5:1 (200 ug:4*10^13)	Yellow, Pink

- 1. Calculate volume of SWNTs needed (stock=20 ug/ml)
- 2. Mix in dialysis tubes (label clips of your tubes)
- 3. Dialyze against NaCl pH 5.3 then 10

Low pH = minimize electrostatic repulsion (phage/SWNT High pH = stabilize complex, ready for TiO<sub>2</sub>

### Guidelines for your 20.109 research proposal

Writing a research proposal requires that you identify an interesting topic, spend lots of time learning about it, and then design some clever experiments to advance the field. It also requires that you articulate your ideas so any reader is convinced of your expertise, your creativity and the significance of your findings, should you have the opportunity to carry out the experiments you've proposed.

### FNT:

- 5 recent findings
- 3-5 sentence description of each
- reference